INSTRUCTION MANUAL

**LCD Professional Digital Microscope
Model # 44345**

Introduction

Thank you for purchasing the Celestron LCD Digital Microscope with a 3.5" touch screen monitor. Your microscope is a precision optical instrument, made of the highest quality materials to ensure durability and long life. It is designed to give you a lifetime of pleasure with a minimal amount of maintenance. This microscope was designed and intended for those age 13 and above.

Before attempting to use your microscope, please read through the instructions to familiarize yourself with the functions and operations to maximize your enjoyment and usage. See the microscope diagram to locate the parts discussed in this manual.

This microscope provides high powers from 40x up to 400x (up to 1600x with digital zoom). This microscope is mainly suited for examining specimen slides of yeasts and molds, cultures, plant and animal parts, fibers, bacteria, etc. You can also examine thin, small objects such as coins, stamps, PC boards, insects, and other objects especially at the lower powers but remember the lowest power is 40x.

The LCD Digital Microscope does not use eyepieces that are used in traditional microscopes. You will view specimens or objects on the LCD screen which are easy to see and you can enjoy them with others also. You can also take snapshots or short videos with the built-in digital camera.

The final section provides simple care and maintenance tips for you to follow to ensure that your microscope provides you with years of quality performance, usage, and enjoyment as well as some troubleshooting hints if you have problems seeing an image.

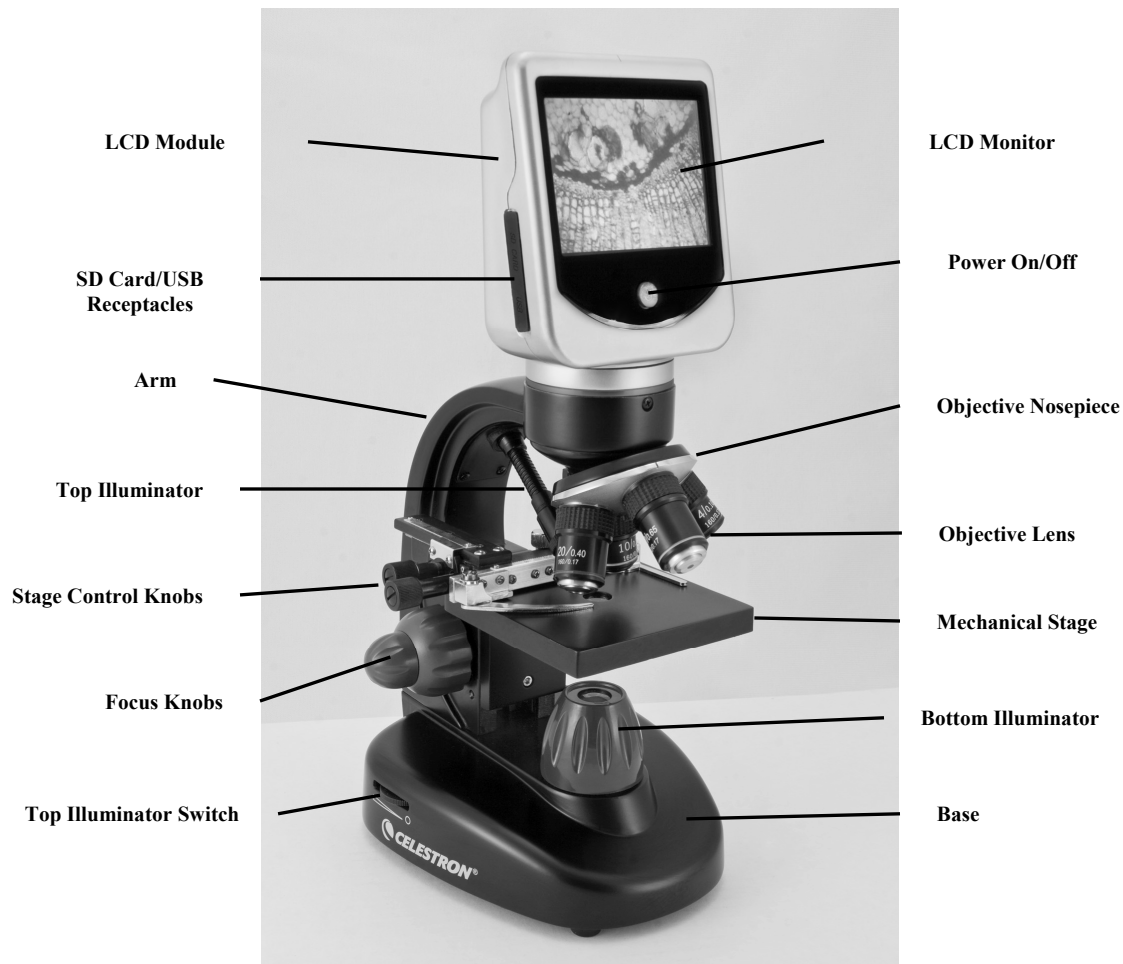
Figure 1

Standard Accessories Included with your Microscope

• Digital Camera – built-in	• Filter Wheel/Diaphragm – 6 position
• LCD Touch Screen Monitor	• AC Adapter
• 4x Objective Lens	• USB Cable 2.0
• 10x Objective Lens	• 7 pcs Prepared Slides
• 20x Objective Lens	• SD Card Slot
• 40x Objective Lens	• Dust Cover
• Top & Bottom Illuminators	• Soft Case

Specifications

Stage	Mechanical Stage 3.5" x 3.5" (88mm x 88mm)
Digital Camera	2MP CMOS; 10x Magnification in lieu of an Eyepiece; 1600x 1200 pixel array
LCD Monitor	3.5" (90mm) with 4x Digital Zoom – High definition touch digital TFT display Resolution – 320 x 240 pixels, Contrast ratio – 300, Color Support – 262K colors, Brightness/Luminance – 430cd/m2
Focuser	Dual – Coarse/Fine
Objectives	Achromatic 4x, 10x 20x, and 40x You can use optional objective lenses up to 60x but not lower than 4x – DIN35 RMS thread size WJ 4/5" x 1/36"
Flash Memory	1GB (approximately 1,300+ snapshots @ 2mp or video of 4 hours+) Snapshots in JPEG format and 3GP files for Video
LCD Rotation	180° -- 90° left / 90° right
Filter Wheel	Red/Green/Blue/1mm hole/3mm hole/ 6mm hole (default)
Nosepiece	Quad with click stop
Illuminators	Built-in electric – both are LED 6Volt and 6Watt
Condenser	N.A. 0.65
AC Adapter	Input Universal 100 to 240Volt 50/60HZ
Batteries	User supplied 4AA – approx. 3 hour usage
Weight/Dimensions	60oz/1.7Kg 6.9" (175mm) x 5.4" (137mm) x 12.7" (323mm)

Magnification (Power) Table

Use the following table to determine the magnification of the different objective lenses in combination with your microscope using the normal mode of the digital image on the LCD screen and using the digital zoom feature.

Objective Lens	4x	10x	20x	40x
Digital Image -- normal	40x	100x	200x	400x
Maximum with 4x digital zoom feature	160x	400x	800x	1600x

Setting Up Your Microscope

1. Carefully remove the microscope and other parts from the carton and set them on a table, desk, or other flat surface.
2. Remove the plastic cover from the microscope.
3. Plug the small cable from the AC adapter into the socket on the back of the base (see Figure 2).
4. Insert the plug you need (see note below) into the AC adapter and then plug the adapter into the proper power source.

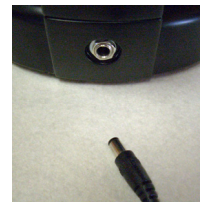
Figure 2

Figure 3

Note: The AC Adapter supplied with this microscope has an interchangeable plug system that can work in the USA, UK, Europe, Australia, and many other parts of the world. You can easily change plugs for your particular requirement by doing the following:

- A. Pull the plug up and away while holding the base of the AC Adapter and the plug will come off.
- B. Take the plug style that you want to use and align the two holes in the bottom of the plug with the two pins in the AC Adapter (see Figure 3). Then, push the plug into the AC Adapter all the way and you are finished.

Battery Operation – You can use your microscope without AC power if you choose. This gives you the freedom to operate the microscope outdoors or indoors anywhere you please. The battery operation requires 4AA batteries (user supplied). Open the battery door on the bottom of the microscope and insert the batteries according to the battery polarity shown in the battery compartment (Figure 4). After the batteries are installed close the battery door. Battery life will typically be three hours.

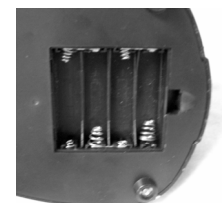
Figure 4

Microscope Operation

Before looking at specimens you must turn the LCD on, turn on the proper illumination, and understand how to use the mechanical stage and then you are ready to begin viewing.

LCD Module

This digital microscope is different than traditional microscopes --- instead of using eyepieces to look at a specimen in a traditional microscope, the LCD monitor replaces the eyepieces so you can look at the specimen on the screen by yourself or share the views with others. To begin to view specimens with your microscope, you will have to turn the LCD monitor on by pushing and holding the Power Button (see Figure 1) until the LED comes on below the on/off button and you see "Celestron Digital Microscope" on the screen. That is basically all you need to do to use the LCD screen for viewing specimens. The touch screen functions on the LCD Module are mainly used for taking images (snapshots and video) and performing other functions and will be discussed later in this manual.

Illumination

To get the sharpest and best views, the proper illumination (lighting) must be chosen:

1. To turn the illuminator(s) on, see Figures 5 & 6 and turn the switches as shown for each.
2. The top illuminator (Figure 1) was designed to be used at low power (4x objective) as higher power objective lenses (10x, 20x & 40x) will block some of the light. If you need to use high power to observe solid objects, use a bright secondary light (desk lamp, etc.) for directed illumination.
3. The bottom illuminator (Figure 1) is used mainly for specimen slides where the light shines up through the hole in the stage through the slide.
4. Having both illuminators on at the same time can provide enough light for thick and irregular specimens.

Figure 5

Adjusting the Lighting

Figure 6

Specimens of different size, thickness, and color variations will require different levels of illumination. Adjust the brightness by turning the switches shown in Figure 5 & 6.

When viewing a specimen that is not transparent or dark in color, you may need to increase the amount of light to resolve certain features or details. This is best done by simply increasing the brightness of the illuminator by rotating the brightness control dial all the way to its highest setting.

Optimum lighting will be found by experimenting with adjustments as each specimen may require slightly different illumination as well as the same specimens viewed under different powers.

Viewing a Specimen

Your instrument is provided with a mechanical stage with a stage holder clamp and directional knobs –see Figure 7 below.

1. Use the clamp lever to open the clamping arm of the stage holder clamp.
2. Place a specimen slide (1" x 3"/25.4mm x 76.2mm size) inside the holder and gently close the clamping arm against the slide.
3. Use the stage movement knobs to position the specimen over the opening in the stage. The rear stage movement knob moves the X axis (forward and backward) whereas the front stage movement knob moves the Y axis (side to side).

Note: A vernier scale on both axes allows the exact marking and replication of an object in the field of view that the user may want to come back to.

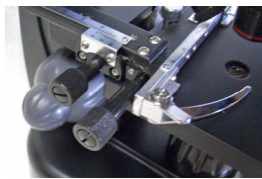
Figure 7

4. Use the objective nosepiece (Figure 1) to rotate the objective lenses (Figure 1) until the 4x objective lens is directly over the specimen. Always start with the lowest power objective (4x with this microscope) which gives you 40 power and work your way up to higher powers. At 40 power you will have the widest field of view and the brightest image.
5. Look at the LCD screen while turning the focus knob (Figures 1 & 7) until the specimen comes into view. You may need to adjust the stage movement knobs (Figure 7) slightly to center the specimen in the field of view. The larger focus knob is the coarse focus and the smaller knob is for fine (exact) focusing.

6. With the 4x objective lens, you can also vary the power anywhere from 40x to 160x by using the digital zoom.
7. For higher powers, you will need to rotate the objective nosepiece to the 10x or 20x and to the 40x objective for the maximum power. You will have to refocus when changing the power of the objective lenses. While using any of these objective lenses you also can increase power by using the digital zoom. Note that using a higher power objective lens will yield sharper images versus a lower power objective lens and digital zoom for the same magnification.
8. To use the **digital zoom**, you touch the screen and push lightly on the right side to increase the power from 1x to 4x or touch the screen and push lightly on the left side to decrease the power.

Note: When changing objective lenses, lower the stage to its lowest position so you will not hit anything during the rotation. Also, at the higher powers, be careful when raising the stage close to the objective lens so that the objective does not hit the slide specimen (or other object) and cause damage.

Using Filters & the Diaphragm

Normally most viewing or imaging will be done without colored filters and before using the microscope check to make sure no filters are in the optical path. However, to bring out different levels of detail, experiment with changing the color of the back lighting of the specimen especially for very bright transparent specimens. To change the lighting color, rotate the wheel (Figure 8 on the right side) to the desired color – Red (R), Green (G), Blue (B). Each color is centered when you hear/feel the faint click stop. You may need to refocus by adjusting the focus knob (Figure 1) slightly for best viewing. You should experiment with each of the colors to see the results.

Diaphragm – within the wheel are holes with three different diameters, (1) 1mm, (3) 3mm, (6) 6mm which limit the amount of light passing through to the specimen. These holes are part of the diaphragm which allows you to change the hole opening size to maximize the contrast, brightness, etc.

The default setting is the (6) for the 6mm hole which should be used for most viewing. You can look under the stage to make sure the proper setting you wish is actually being used.

Wheel

Figure 8

Figure 9

Rotating the LCD Screen

You can rotate the viewing position of the LCD screen 180° -- 90° to the right and 90° to the left. You can view any position you choose along the 180° rotation. This function allows you to share the view with others without actually moving the complete microscope. To move the LCD screen, hold the round black part (see Figure 1) below the orange ring with one hand and then hold the LCD module with the other and move it to the position you desire.

Turn the LCD screen on and you are now ready to use your microscope for viewing and if any problems check the trouble shooting section.

Digital Imaging

You can take snapshots or a short video with your microscope with the built-in digital camera. With the internal storage memory, you do not need to use a PC or any other devices to do imaging. Transferring of images to your PC for saving and or printing them is easy and will be discussed later in this manual.

Note: If you are going to take images, do not connect the USB cable to your PC or damage could occur to the equipment. The USB cable is not used at all for taking images.

Settings and Information for the Digital Camera and the Touch Screen

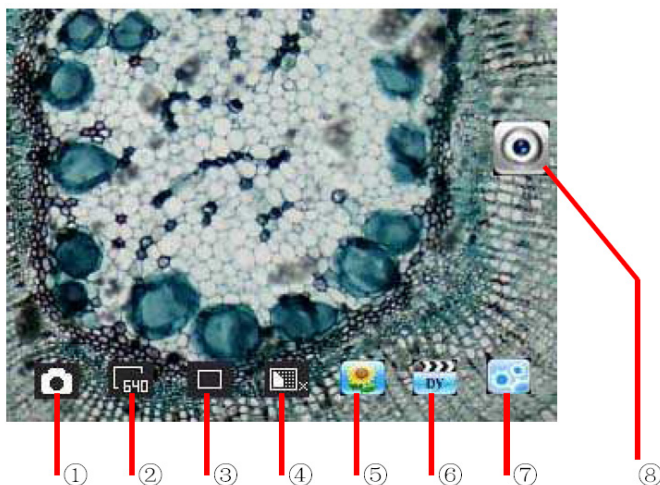
Figure 10

The touch screen icons and their functions are quite easy to use and intuitive in nature. Below will be discussed the general use of the icons.

The touch screen has various functions and choices among those functions. From the image to the left you will find the following eight icons:

1. Video mode
2. Pixel setting
3. Snapshot selection
4. Color effect
5. Image review
6. Video review
7. Settings
8. Viewing / Snapshot mode

Note: After turning the LCD Monitor off, most settings will revert to the default settings

Setting icon (7)

Language – choose from Chinese (simple or traditional), English, French, German, Italian, Japanese, Korean, Portuguese, Russian, and Spanish

Time for Auto Play Review – 1 second to 6 seconds

Memory – internal memory or SD Card

Time & Date – year, month, and date

Calibration – touch screen sensitivity

Brightness – of touch screen

Auto Shutdown of Power – never to 5 minutes

Default Factory Settings – change all back to factory settings

Pixel setting icon (2)

Lightly tap the icon to change the pixel setting from 320 (320x640), 640 (640x480), 1280 (1280x1024), 1600 (1600x1200), 2048 (2048x1536), and 2560 (2560x 2048). 1600 is the sensor resolution and 2048 and 2560 give you higher resolution through interpolation.

Snapshot selection (3)

Normal is single shot but you can set for timed or a burst of three shots.

Color effect (4)

Lightly tap the icon to change from black & white, sepia, negative, transparent, and normal (x).

Viewing / Snapshot mode (8)

When viewing specimens, you will see this icon on the right of the screen. After you have made the various settings above and you want to take snapshot images, you just lightly press the icon to take an image.

Image review (5)

In this mode, you can review the snapshot images you have taken. Just touch the arrows and touch the screen to navigate and to see full screen images, etc.

Video mode (1)

You can take videos in this mode.

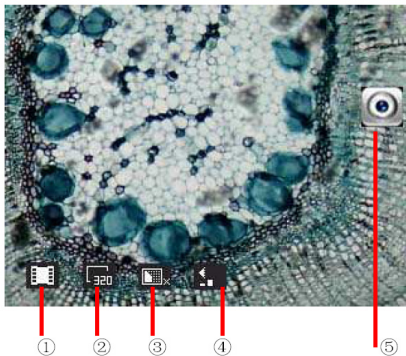
Figure 11

To take videos, you need to make the settings first. In the image to the left, you will see the following icons on the screen after you press the Video icon. The icons on this image (Figure 11) do the following:

1. Touch to go back to the viewing mode
2. Set the pixel size to 160 or 320 (320 being a higher resolution for videos) by touching the screen
3. Set the color desired – black & white, sepia, negative, or normal by touching the screen
4. Video quality setting – low, medium, or high quality by touching the screen
5. Record Video – touch to begin video and touch again to stop the video.

Video review (6)

In this mode you can review the videos you have taken. Just touch the arrows and touch/scroll the screen to navigate and view your videos and delete them if you choose.

Using an SD Card

You can also take images (snapshot or video) by using an SD (Secure Digital) Card (4GB max). Your SD card is inserted in the LCD Monitor (Figure 1).

When you insert the SD card properly the SD card icon will be visible on the LCD screen. When imaging this way the images are saved on the SD card and not in the internal storage memory.

Note: Inserting or removing an SD card while the LCD is on may cause the LCD to shut down and/or could damage the SD card.

Transferring Your Images

To transfer images to a PC or MAC, you need to have a free USB port and have an imaging program for snapshots and/or video.

Note: Do not disconnect the USB cable while transferring images or damage may occur.

1. You can transfer images from the internal storage memory to your PC by using the supplied USB Cable. The small plug end of the cable plugs into the LCD Monitor (see Figure 1)) and the large plug end of the cable plugs into your PC. If the connections are proper you will see on your microscope screen “MSDC” or similar data. Your PC will automatically recognize the new hardware. Then you will choose which program on your PC you want to transfer the images to.
2. If you used an SD card to store your images, you can transfer them to your PC using the method in # 1 above or you can take the SD card out and use the SD card slot on your PC to transfer the images. In either way your PC will ask you to choose which program you want to transfer the images to.

Trouble Shooting

If you do not get an image to view on your LCD screen, here are a few things to double check:

1. Make sure the AC Adapter is plugged in to an AC power source and attached to the microscope securely and correctly.
2. Make sure you have the illuminator turned on with maximum brightness adjustment.
3. Make sure the objective lens you have chosen is set correctly and it has clicked in the right position.
4. Make sure that the diaphragm (filter wheel) is set correctly at a click position so that the illuminated light comes up properly – the normal position is the 6 position (6mm hole) for most applications.
5. Make sure the specimen slide is correctly fit into the clamp on the mechanical stage and properly centered.

Care, Maintenance, and Warranty

Your Celestron microscope is a precision optical instrument and should be treated with care at all times. Follow these care and maintenance suggestions and your microscope will need very little maintenance throughout its lifetime.

- When you are done using your microscope, remove any specimens left on the stage.
- Turn off the illuminator switches.
- Turn off the LCD monitor – push and hold the on/off button until you see “Power Shutting Off”.
- Unplug the power cord.
- Always place the plastic bag or dust cover over the microscope when not in use or when being stored to help keep it clean.
- Store the microscope in a dry and clean place.
- Be very careful if using your microscope in direct sun light to prevent damage to the microscope or your eyes.
- When moving your microscope, carry it by the “arm” with one hand and not by the focuser knob, LCD monitor, etc. Then, put your other hand under the base for support.
- Clean the outside surfaces (metal and plastics) with a moist cloth.
- Always unplug any cords before cleaning.
- Never clean optical surfaces with cloth or paper towels as they can scratch optical surfaces easily.
- Blow off dust with a camel’s hair brush or an air blower from optical surfaces.
- To clean fingerprints off of optical surfaces, use a lens cleaning agent and lens tissue available at most photo outlets and when cleaning do not rub in circles as this may cause streaks or scratches to occur.
- Never disassemble or clean internal optical surfaces. This should be done by qualified technicians at the factory or other authorized repair facilities.
- When handling glass specimen slides, use care as the edges can be sharp.

Warranty

Your microscope has a two year limited warranty. Please see the Celestron website for detailed information at www.celestron.com.

EEC: This product complies with EEC guidelines in EN61558-2-6:1997 and EN61558-1:1997+A1

FCC Statement

This device complies with Part 15 of FCC Rules. Operation is subject to the following two conditions:

1. This device may not cause harmful interference, and
2. This device must accept any interference received, including interference that may cause undesired operation.

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Torrance, California 90503 U.S.A.